

Luria Broth (Miller's)

Cat. # P1310-1000, P1310-5000

Also Known as:	N/A
Form:	Granular
Quantity:	See tube label
Technical Specification:	Components in 25 g mixtures: Peptone: 10 g Yeast Extract: 5 g Sodium Chloride: 10 g
Description:	Mixtures of peptone, yeast extract and NaCl that support bacterial growth. Compared to powder, granulated formula can significantly reduce dust during weighing in preparation of culture.
Storage:	Ship under ambient. Keep container tightly closed and store at a dry place under room temperature.
Usage	Instruction for preparing Luria Broth culture: Dissolve 25 g in 1 L distilled water in a glass flask, heat and agitate to completely dissolve, adjust pH to 7.0, then autoclave at 121 °C for 15 min.